

TACHOM-UN

Method of Use

The QLA Handheld Point-And-Shoot Tachometer (TACHOM-UN) targets the rotating Basket or Paddle shaft and provides instant precision results.

Measurements are made by placing a small cutting of reflective tape (approx 1/8" X 1/2" Long) lengthwise on the rotating basket or paddle shaft (Supplied with purchase of TACHOM-UN). Point the Tachometer at the strip of reflective tape. The Red LED light locks on the self-adhesive reflective tape and counts the rotations of the shaft. Within two seconds it provides your reading from 10 to 99,999 revolutions per minute (RPM).

